GalactEXO[™]

FOR RESEARCH USE ONLY STORE AT



SmartEnzymes[™]





INSTRUCTIONS FOR PRODUCT

GalactEXO 2000 units (G1-GM1-020) Digestion of up to 2 mg glycoprotein

Last revised July 2020

QUICK GUIDE



 Reconstitute GalactEXO in 100 µl ddH₂O to a concentration of 20 units/µl



 Add 1 unit GalactEXO / 1 µg glycoprotein or 1 pmol oligosaccharides



Incubate for 2 h to 18 h at 37°C



GalactEXO is a mix of β -galactosidases for efficient removal of galactose residues on Nand O-glycosylated proteins or oligosaccharides. The mix is composed of two galactosidases for highly efficient hydrolysis of both β 1-3 and β 1-4 linked galactoses¹. GalactEXO hydrolyzes glycoproteins under native conditions and displays a high activity in a broad pH range, 5.5 to 7.5. The enzymes in GalactEXO are derived from *Akkermansia muciniphila* and expressed in *E. coli*. The two galactosidases have his tags and the molecular weights of the components are 87 kDa and 109 kDa, respectively.

Unit Definition

One unit of GalactEXO hydrolyzes galactoses from >95 % of 0.15 nmol Gal- β 1,3-GalNAc-pNP and 2 nmol Gal- β 1,4-GlcNAc-pNP when incubated in 20 mM Tris pH 6.8 at 37°C for 30 min.

Content and Storage

GalactEXO is supplied lyophilized in TBS pH 7.6, with no preservatives added.

GalactEXO is shipped cold and should be stored at -20°C upon arrival.

After reconstitution GalactEXO is stable for 1 month at $+4-8^{\circ}$ C.

GalactEXO is for R&D use only.

DETAILED PROTOCOL

Additional Materials Required

• Reaction buffer²: 20 mM Tris pH 6.8

Sample Preparation

• Prepare the glycoprotein of interest in the reaction buffer at a concentration of 0.5-5.0 mg/ml.

Digestion of Glycoproteins

1 Prepare GalactEXO

 Reconstitute GalactEXO in 100 µl ddH₂O to a concentration of 20 units/µl

2 Add GalactEXO

Add 1 unit GalactEXO / 1 µg glycoprotein³

3 Digestion

Incubate for 2 to 18h at 37°C

Optimization of enzyme concentrations and incubation time may be needed depending on the substrate.



Digestion of Oligosaccharides

1 Prepare GalactEXO

 Reconstitute GalactEXO in 100 µl ddH₂O to a concentration of 20 units/µl

2 Add GalactEXO

 Add 1 unit GalactEXO / 1 pmol oligosacharides

3 Digestion

Incubate for 1 h at 37°C

Notes

1. GalactEXO also hydrolyses β1-6 linked galactoses to a certain extent.

 GalactEXO displays high activity in buffers with pH values from 5.5 to 7.5 and over a wide range of ionic strength. Some optimizations might be required if a buffer other than the recommended reaction buffer is used.

A higher enzyme concentration may increase digestion efficiency of individual glycoproteins. This requires optimization.

Quality Control

GalactEXO is tested to meet the specifications and lot-to-lot consistency.

GalactEXO is tested for the absence of microbial contamination using blood agar plates, Sabouraud dextrose agar plates and fluid thioglycollate medium.

Related Products

SialEXO[®]

For complete removal of α 2-3, α 2-6 and α 2-8 linked sialic acids.

SialEXO[®] 23

For complete and specific removal of α2-3 linked sialic acids.

Immobilized SialEXO®

Complete removal of sialic acids in a spin column format.

Legal and Disclaimers

All rights reserved. Genovis products may be covered by one or more patents, trademarks and copyrights owned or controlled by Genovis AB.

For more information about commercial rights, please contact Genovis team at info@genovis.com

Genovis products are intended for research use only. They are not intended to be used for therapeutic or diagnostic purposes in humans or animals.

© Genovis 2020

SialEXO°

Efficient removal of sialic acids on O-glycosylated and N-glycosylated proteins

SialEXO is a mix of sialidases for efficient removal of sialic acids on O-glycosylated and N-glyco-sylated proteins. The mix is composed of two sialidases for highly efficient hydrolysis of α 2-3, α 2-6 or α 2-8 bonds.

SialEXO hydrolyzes glycoproteins under native conditions and displays a high activity in a broad pH range, 6.5 to 9.

The enzymes in SialEXO are derived from Akkermansia muciniphila and expressed in E. coli. SialEXO is composed of two sialidases with His-tags and the molecular weights of the components are 42.8 kDa and 65.7 kDa, respectively.



SialEXO[®] 23

Specific removal of α2-3 linked sialic acids on O- and N-glycans

SialEXO 23 is a sialidase for specific removal of α 2-3 linked sialic acids on O- and N-glycans.

SialEXO 23 hydrolyzes sialic acids on glycans under native conditions and displays a high activity in a broad pH range, 7 to 9. The enzyme is derived from Akkermansia muciniphila and expressed in E. coli. The enzyme contains a His-tag and the molecular weight is 66 kDa.

SialEXO 23 can be used for exoglycosidase sequencing by combining results from a reaction with the sialidase product SialEXO, that efficiently hydrolyzes α 2-3, α 2-6 and α 2-8 linked glycosidic bonds.





USA & Canada

Genovis Inc. 245 First Street, Suite 1800 Cambridge, MA 02142 USA

Customer service: 1-617-444-8421 Order phone (toll free): 1-855-782-0084 Order fax: 1-858-524-3006 Email: orders.us@genovis.com

EMEA & Asia

Genovis AB Box 790 SE-220 07 Lund Sweden

Customer service: 0046 (0)46 10 12 30 Order phone: 0046 (0)46 10 12 30 Order fax: 0046 (0)46 12 80 20 Email: order@genovis.com