

GlySERIAS™

FOR RESEARCH
USE ONLY
www.genovis.com

STORE AT
-20°C



SmartEnzymes™



GENOVIS

INSTRUCTIONS FOR PRODUCT

GlySERIAS, 2000 units (A0-GS1-020)

Digestion of up to 2 mg fusion protein

Quick Guide

1 Prepare GlySERIAS™

- Reconstitute GlySERIAS in 50 μ l ddH₂O to a concentration of 40 units/ μ l.




2 Add GlySERIAS™

- Add 1 unit GlySERIAS / 1 μ g fusion protein.



3 Digestion

- Incubate for 1 h at 37°C.

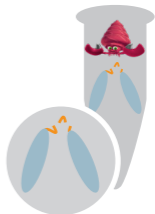

PRODUCT DESCRIPTION

GlySERIAS is an enzyme that digests flexible linkers in engineered fusion proteins containing two or more proteins or peptides. It is active on $(\text{Gly}_4\text{Ser})_n$, Gly_xSer_y and polyglycine linkers. The repetitive design of the linker will lead to several simultaneous digestion sites and separation of the previously linked components. Optimal activity occurs at 37°C, pH 7.6 under native conditions.

GlySERIAS is cloned from phage K and is recombinantly expressed in *E. coli*. The enzyme contains a His-tag and has a molecular weight of 18 kDa.

Unit Definition

One unit GlySERIAS digests $\geq 95\%$ of 1 μg dulaglutide at a minimum of one site, when incubated in TBS (50 mM Tris-HCl, 150 mM NaCl, pH 7.6) at 37°C for 15 minutes.

Content and Storage

GlySERIAS is supplied lyophilized in 50 mM Tris-HCl, 150 mM NaCl, pH 7.6, with no preservatives added. Upon arrival, the enzyme should be stored at -20°C.

After reconstitution, the GlySERIAS enzyme is stable for at least 1 month at +4-8°C.

GlySERIAS is for R&D use only.

DETAILED PROTOCOL

Additional Materials Required

- Digestion buffer: TBS, pH 7.6.

Sample Preparation

Prepare the fusion protein in the digestion buffer¹. The final protein concentration in the digestion reaction should be 1-5 mg/ml.

1 Prepare GlySERIAS™

- Reconstitute GlySERIAS in 50 µl ddH₂O to 40 units/µl.

2 Add GlySERIAS™

- Add 1 unit GlySERIAS / 1 µg fusion protein².

3 Digestion

- Incubate for 1 h at 37°C^{3,4}.

Notes

- 1. Optimal activity is achieved in TBS buffer pH 7.6. The enzyme is active in pH 6.5-9 but the digestion efficiency may differ between different GS-linked proteins.*
- 2. A higher enzyme concentration may increase digestion efficiency of individual GS-linked proteins. This requires optimization.*
- 3. A shorter incubation time will allow for a more complete coverage of linker sequence whereas a longer incubation time will reduce complexity and result in more homogeneous subunits.*
- 4. The linker may not be completely removed from the GS-linked proteins.*

Quality Control

GlySERIAS is tested to meet the specifications and lot-to-lot consistency.

GlySERIAS is tested for absence of microbial contamination with blood agar plates, Sabouraud dextrose agar plates and fluid thioglycollate medium.

Related Products

FabRICATOR®

Below hinge digestion of IgG

FabALACTICA®

Above hinge digestion of human IgG1

FabDELLO™

Above hinge digestion of human IgG1

Legal and Disclaimers

All rights reserved. Genovis products may be covered by one or more patents, trademarks and copyrights owned or controlled by Genovis AB. For more information about commercial rights, please contact the Genovis team at licensing@genovis.com.

Genovis products are intended for research use only. They are not intended to be used for therapeutic or diagnostic purposes in humans or animals.

All goods and services are sold subject to Genovis' General Terms and Conditions of Sale.

© Genovis AB.



USA & Canada

Genovis Inc.
245 First Street, Suite 1800
Cambridge, MA 02142
USA

Customer service: 1-617-444-8421
Order phone (toll free): 1-855-782-0084
Order fax: 1-858-524-3006
Email: orders.us@genovis.com

EMEA & Asia

Genovis AB
Box 790
SE-220 07 Lund
Sweden

Customer service: +46 46 10 12 30
Order phone: +46 46 10 12 30
Order fax: +46 46 12 80 20
Email: order@genovis.com